

## Integrity Test Unit BECO® Control Mini



The BECO Control Mini integrity test unit from Eaton's Begerow Product Line facilitates the execution of integrity tests. Because the measurement of the pressure drop during the test period is extremely precise, the exact diffusion rate in ml/min can be determined. At the end of the integrity test, the measuring instrument indicates whether the measurement result was positive or negative. Every test is recorded and may be viewed over the display or via data cable over PC/laptop (Version < Windows 7) so that results can be documented.

By entering individual filter data, any cartridge filter or other filter media can be evaluated using this measuring instrument. For this purpose, 19 programmable test programs are available.

Values to be entered, one-time input:

- Pore size
- Diffusion rate/cartridge allowed
- Test pressure
- Number of cartridges
- Filter volume
- Stabilization
- Duration of test

### Benefits

- Accurate, reproducible results
- Portable application
- Easy handling
- Reliable documentation

### Accomplishments

- High accuracy of the pressure sensor and permanent error diagnostic
- Menu-driven prompt
- High memory capacity and user-friendly data transfer
- Encapsulated and splash waterproof design (IP 64)



BECO Control Mini with interface unit

### Accessories for BECO Control Mini

Article number	Description
BJTG1D20	Control Mini
BJTG1025	Interface unit Control Mini
BJTG3D26	Rectus-coupling with sleeve nut

#### **North America - HQ**

44 Apple Street,  
Tinton Falls, NJ 07724  
Toll Free: 800 656-3344  
(North America Only)  
Voice: +1 732 212-4700

#### **Europe/Africa/Middle East**

Auf der Heide 2  
53947 Nettersheim, Germany  
Voice: +49 2486 809-0

#### **Internormen Product Line**

Friedensstraße 41  
68804 Altlusheim, Germany  
Voice: +49 6205 2094-0

#### **Begerow Product Line**

An den Nahewiesen 24  
55450 Langenlonsheim, Germany  
Voice: +49 6704 204-0

#### **Brazil**

Av. Julia Gaioli, 474 - Bonsucesso  
07251-500 - Guarulhos  
Brazil  
Voice: +55 11 2465 8822

#### **China**

No. 7 Lane 280 Linhong Road,  
Changning District,  
Shanghai 200335, China  
Voice: +86 21 5200 0422

#### **Singapore**

4 Loyang Lane #04-01/02  
Singapore 508914  
Voice: +65 6825 1668

For more information, please e-mail us at [filtration@eaton.com](mailto:filtration@eaton.com) or visit us online at [eaton.com/filtration](http://eaton.com/filtration) for a complete list of Eaton's filtration products.

Not all products in Eaton's Begerow Product Line are available in all regions. Please contact your local Eaton office to determine availability.

© 2012 Eaton Corporation. All Rights Reserved. All trademarks and registered trademarks are the property of their respective owners.

All information and recommendations appearing in this brochure concerning the use of products described herein are based on tests believed to be reliable. However, it is the user's responsibility to determine the suitability for his own use of such products. Since the actual use by others is beyond our control, no guarantee, expressed or implied, is made by Eaton as to the effects of such use or the results to be obtained. Eaton assumes no liability arising out of the use by others of such products. Nor is the information herein to be construed as absolutely complete, since additional information may be necessary or desirable when particular or exceptional conditions or circumstances exist or because of applicable laws or government regulations.



*Powering Business Worldwide*