

Depth Filtration
BECO® PR Range



Depth Filter Sheets for the Pharmaceutical Industry

BECO PR depth filter sheets from Eaton's Begerow Product Line meet the high demands of the pharmaceutical industry. Exceptionally pure raw materials and a special production method produce BECO PR depth filter sheets with low endotoxin content. The special characteristic of this range is high endotoxin retention during the filtration of many pharmaceutical products.

The specific advantages of BECO PR depth filter sheets:

- High endotoxin retention as well as a maximum microbe retention rate.
- The innovative production process guarantees an endotoxin content of less than < 0.125 EU/ml.
- Maximum raw material purity for minimum migration of soluble ions.
- The ideal combination of various filtration mechanisms (surface, adsorption, depth filtration) and adsorptive properties ensures maximum reliability.
- Comprehensive quality assurance for all raw and auxiliary materials and intensive in-process controls ensure consistent quality of the finished products.
- Before delivery, the endotoxin content of < 0.125 EU/ml of all BECO PR depth filter sheets is tested with the help of an LAL test. A certificate is available on request.

Microbe Reduction and Removal

BECO PR Steril S 100, PR Steril S 80, PR Steril 40
BECO depth filter sheets boast high microbe retention rates achieved through the tight-pored structure and an electrokinetic potential with an adsorptive effect.

These depth filter sheets are particularly suitable as prefilters for subsequent membrane filtration because of their high capacity to retain endotoxins and colloidal components.

Fine Filtration


BECO PR 12

BECO depth filter sheets for achieving a high degree of clarification. These depth filter sheets reliably retain ultra-fine particles and provide bioburden reduction.

In practice, these depth filter sheets serve as ideal prefilters for protection of membrane filters, reverse osmosis systems, and to protect chromatography columns.



Water throughput BECO PR range

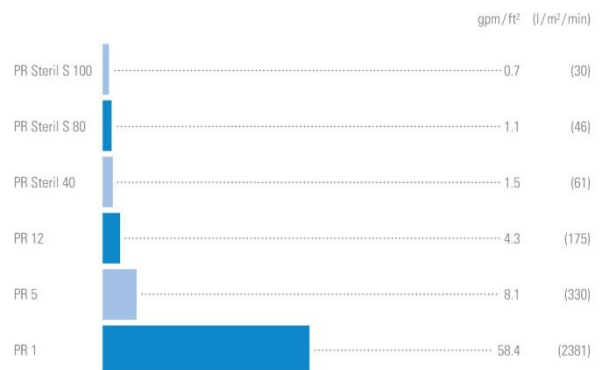

Conditions: Δ p = 14.5 psi (100 kPa, 1 bar), Medium: Water at 68 °F (20 °C)

Clarifying Filtration and Coarse Filtration

BECO PR 5, PR 1

BECO depth filter sheets with large-volume cavity structure. These depth filter sheets have a high dirt holding capacity for particles and are very suitable for clarifying filtration applications.

Physical Data

This information is intended as a guideline for the selection of BECO depth filter sheets.

Type	Article no.	Nominal retention rate µm	Thickness in (mm)	Ash content %	Bursting strength wet		Water throughput at		Endotoxin content** EU/ml
					psi	(kPa)	Δ p = 14.5 psi gpm/ft ²	(Δ p = 100 kPa* l/m ² /min)	
PR Steril S100	27295	0.1	0.15 (3.9)	58.0	> 7.3	(50)	0.7	(30)	< 0.125
PR Steril S80	27280	0.2	0.15 (3.9)	50.0	> 11.6	(80)	1.1	(46)	< 0.125
PR Steril 40	27240	0.4	0.15 (3.9)	49.0	> 7.3	(50)	1.5	(61)	< 0.125
PR 12	27212	0.8	0.15 (3.9)	50.0	> 18.9	(130)	4.3	(175)	< 0.125
PR 5	27205	2.0	0.15 (3.9)	50.0	> 8.7	(60)	8.1	(330)	< 0.125
PR 1	27200	4.0	0.17 (4.3)	48.0	> 6.5	(45)	58.4	(2381)	< 0.125

The water flow is a laboratory value characterizing the different BECO depth filter sheet types. It is not the recommended flow rate.

* 100 kPa = 1 bar

** Endotoxin content analysis after rinsing with 13.2 gal (50 l) of endotoxin-free water per square meter

Chemical Data

BECO depth filter sheets meet the requirements of LFGB*, Recommendation XXXVI/1 issued by BfR**, and the test criteria of FDA*** Directive CFR 21 § 177.2260.

Chemical resistance of the BECO depth filter sheets to different solvents over a contact time of 3 hours at 68 °F (20 °C).

Solvent	Mechanical strength	Solvent appearance	Solvent	Mechanical strength	Solvent appearance	Solvent	Mechanical strength	Solvent appearance
Aqueous solutions:						Organic solvents:		
Sugar solution, 10%	r	nc	Hydrochloric acid, 1%	r	nc	Methanol	r	nc
With 1% free chlorine	r	nc	Hydrochloric acid, 3%	r	nc	Ethanol	r	nc
With 1% hydrogen peroxide	r	nc	Hydrochloric acid, 5%	r	nc	Isopropanol	r	nc
With 30% formaldehyde	r	nc	Hydrochloric acid, 10%	r	nc	Toluene	r	nc
With 10% ethanol	r	nc	Nitric acid, 1%	r	nc	Xylene	r	nc
With 40% ethanol	r	nc	Nitric acid, 3%	r	nc	Acetone	r	nc
With 98% ethanol	r	nc	Nitric acid, 5%	r	nc	Methyl ethyl ketone	r	nc
Caustic soda, 1%	r	nc	Nitric acid, 10%	r	nc	n-hexane	r	nc
Caustic soda, 2%	r	nc	Sulfuric acid, 1%	r	nc	Dioxan	r	nc
Caustic soda, 4%	r	0	Sulfuric acid, 3%	r	nc	Cyclohexane	r	nc
Ammonia solution, 1%	r	nc	Sulfuric acid, 5%	r	nc	Tetrachloroethylene	r	nc
Ammonia solution, 3%	r	nc	Sulfuric acid, 10%	r	nc	Ethylene glycol	r	nc
Ammonia solution, 5%	r	nc	Acetic acid, 1%	r	nc	Dimethyl sulfide	r	nc
			Acetic acid, 3%	r	nc	N, N-Dimethyl formamide	r	nc
			Acetic acid, 5%	r	nc			
			Acetic acid, 10%	r	0			
<i>r = resistant</i>			<i>nc = no change</i>			<i>0 = slight opalescence</i>		
* = German Food, Commodity, and Feed Act			** = Federal Institute of Risk Assessment			*** = Food and Drug Administration; USA		

Pyrogens/Endotoxins

Pyrogens are biological or chemical substances which can induce a rise in body temperature. One common example are endotoxins. These are cell wall components known as lipopolysaccharides, that are embedded in the outer membrane of gram-negative bacteria.

Quantitative evidence of endotoxins can be determined using the LAL test (**L**imulus **A**mebocyte **L**ysate). This method is an efficient and economical alternative to the rabbit fever test. An independent institute examines the depth filter sheets.

The endotoxin content of the specimens examined is specified in EU/ml (**E**ndotoxin **U**nits). The measurement is carried out after rinsing with 13.2 gal (50 l) of endotoxin-free water per square meter.

Endotoxin Retention Rate

To measure endotoxin retention, a 40% glucose solution containing a defined amount of lipopolysaccharide (LPS) in pyrogen-free water is passed through a depth filter sheet. A defined sample of the filtrate is then measured by means of the LAL test.


Filtration flow rate: 12.3 gal/ft²/h
(500 l/m²/h)

Sampling after: 13.2 gal/sqm and 66 gal/sqm
(50 l/m² and 250 l/m²)

Amount of endotoxin added: 2.2 mg LPS *E. Coli* 055:B5, this equals 4.4 µg LPS/ml or 4.4 x 10⁴ EU/ml

The endotoxin retention rate is indicated in the following graphic.

Endotoxin Retention Rate of BECO PR Depth Filter Sheets

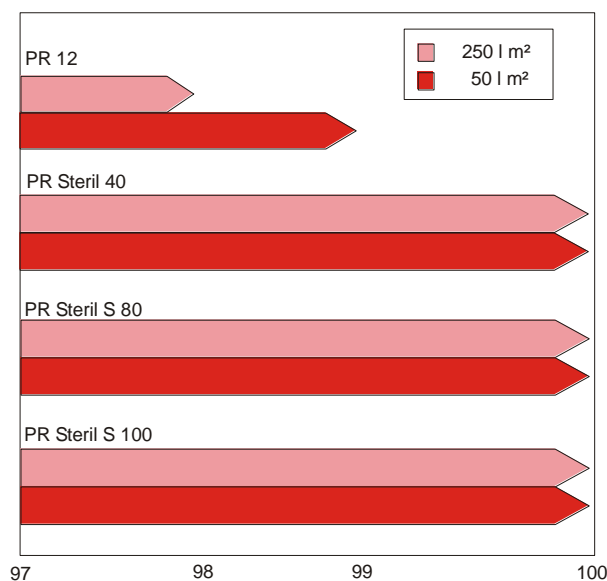

Application Examples:

Application	Germ-reducing filtration	Fine filtration	Clarifying filtration
	PR Steril S 100, PR Steril S 80, PR Steril 40	PR 12	PR 5, PR 1
Dialysis concentrate			x
Human albumin		x	
Photoresist		x	
I-globulin		x	x
Coagulation factors		x	x
Plasma expander solutions	x	x	
Enzyme production		x	x
Hormones	x	x	x
Amino acids	x	x	x
Infusion solutions	x	x	x
Vaccine production	x	x	x
Serums from rabbits, sheep, horses, cattle, calves	x	x	x

Components

BECO depth filter sheets are made from particularly pure natural materials, i.e., finely fibrillated cellulose fibers from deciduous and coniferous trees, cationic charge carriers, and high quality, particularly pure diatomaceous earth.

Instructions for Correct Use

BECO depth filter sheets require careful handling when inserting them into the plate and frame filter. Avoid banging, bending, and rubbing the sheets. Do not use damaged depth filter sheets.

Inserting

Each BECO depth filter sheet has a rough side and a smooth side. The rough side of the depth filter sheet is the unfiltrate side; the smooth side is the filtrate side. Always ensure that the filtrate side is in contact with the clear filtrate plate when inserting the sheets.

Sterilization (Optional)

The wetted BECO depth filter sheets may be sterilized with hot water or saturated steam up to a maximum temperature of **273.2 °F (134 °C)**. The pressed filter package should be loosened slightly. Make sure to sterilize the entire filter system thoroughly. Do not apply final pressure until after the filter package has cooled down.

Sterilizing with Steam

The steam must be free of foreign particles and impurities.

Temperature:	Max. 273.2 °F (134 °C) (saturated steam)
Duration:	Approx. 20 minutes after steam exits from all filter valves.
Rinsing:	After sterilizing with 13.2 gal/sqm (50 l/m ²) at 1.25 times the flow rate.

Filter Preparation and Filtration

Unless already completed after sterilization, Eaton recommends pre-rinsing the closed filter with 13.2 gal (50 l) of water per square meter at 1.25 times the flow rate prior to the first filtration. Depending on the application, this usually equals a rinsing time of 10 to 20 minutes. Test the entire filter for leakage at maximum operating pressure.

High-proof alcohol solutions and chemical products that do not allow pre-rinsing with water should be circulated for 10 to 20 minutes. Dispose of the rinsing solution after rinsing.

Differential Pressure

Terminate the filtration process when a differential pressure of 43.5 psi (300 kPa, 3 bar) is reached. For safety reasons, a differential pressure of 21.8 psi (150 kPa, 1.5 bar) should not be exceeded in applications for separating microorganisms.

Safety

When used and handled correctly, there are no known unfavorable effects associated with this product.

Further safety information can be found in the relevant Material Safety Data Sheet, which can be downloaded from our website.

North America - HQ

44 Apple Street,
Tinton Falls, NJ 07724
Toll Free: 800 656-3344
(North America Only)
Voice: +1 732 212-4700

Europe/Africa/Middle East

Auf der Heide 2
53947 Nettersheim, Germany
Voice: +49 2486 809-0

Internormen Product Line

Friedensstraße 41
68804 Altlussheim, Germany
Voice: +49 6205 2094-0

Begerow Product Line

An den Nahewiesen 24
55450 Langenlonsheim, Germany
Voice: +49 6704 204-0

Brazil

Av. Julia Gaioli, 474 - Bonsucesso
07251-500 - Guarulhos
Brazil
Voice: +55 11 2465 8822

China

No. 7 Lane 280 Linhong Road,
Changning District,
Shanghai 200335, China
Voice: +86 21 5200 0422

Singapore

4 Loyang Lane #04-01/02
Singapore 508914
Voice: +65 6825 1668

Waste Disposal

Due to their composition BECO depth filter sheets are biodegradable. Comply with relevant current regulations, depending on the filtered product.

Storage

BECO depth filter sheets consist of strongly adsorptive materials. The product must be handled carefully during shipping and storage. Store the depth filter sheets in a dry, odor-free, and well-ventilated place.

Do not expose the depth filter sheets to direct sunlight.

BECO depth filter sheets are intended for immediate use and should be used within 24 months of delivery.

Available Formats

All common square or round filter sizes are available for delivery. Special formats are available on request.

HS Customs Tariff: 4812 00 00

Quality Assurance According to DIN EN ISO 9001

Eaton's Begerow Product Line comprehensive Quality Management System has been certified according to DIN EN ISO 9001.

This certification verifies that a fully functioning comprehensive Quality Assurance System covering product development, contract controls, choice of suppliers, receiving inspections, production, final inspection, inventory management, and shipment has been implemented. Extensive quality assurance measures incorporate adherence to technical function criteria and chemical purity and quality recognized as safe under the German legislation governing the production of foods and beverages.

For more information, please e-mail us at filtration@eaton.com or visit us online at eaton.com/filtration for a complete list of Eaton's filtration products.

Not all products in Eaton's Begerow Product Line are available in all regions. Please contact your local Eaton office to determine availability.

© 2012 Eaton Corporation. All Rights Reserved. All trademarks and registered trademarks are the property of their respective owners.

All information and recommendations appearing in this brochure concerning the use of products described herein are based on tests believed to be reliable. However, it is the user's responsibility to determine the suitability for his own use of such products. Since the actual use by others is beyond our control, no guarantee, expressed or implied, is made by Eaton as to the effects of such use or the results to be obtained. Eaton assumes no liability arising out of the use by others of such products. Nor is the information herein to be construed as absolutely complete, since additional information may be necessary or desirable when particular or exceptional conditions or circumstances exist or because of applicable laws or government regulations.



Powering Business Worldwide